


# K series 生物顕微鏡: BIOLOGICAL MICROSCOPES

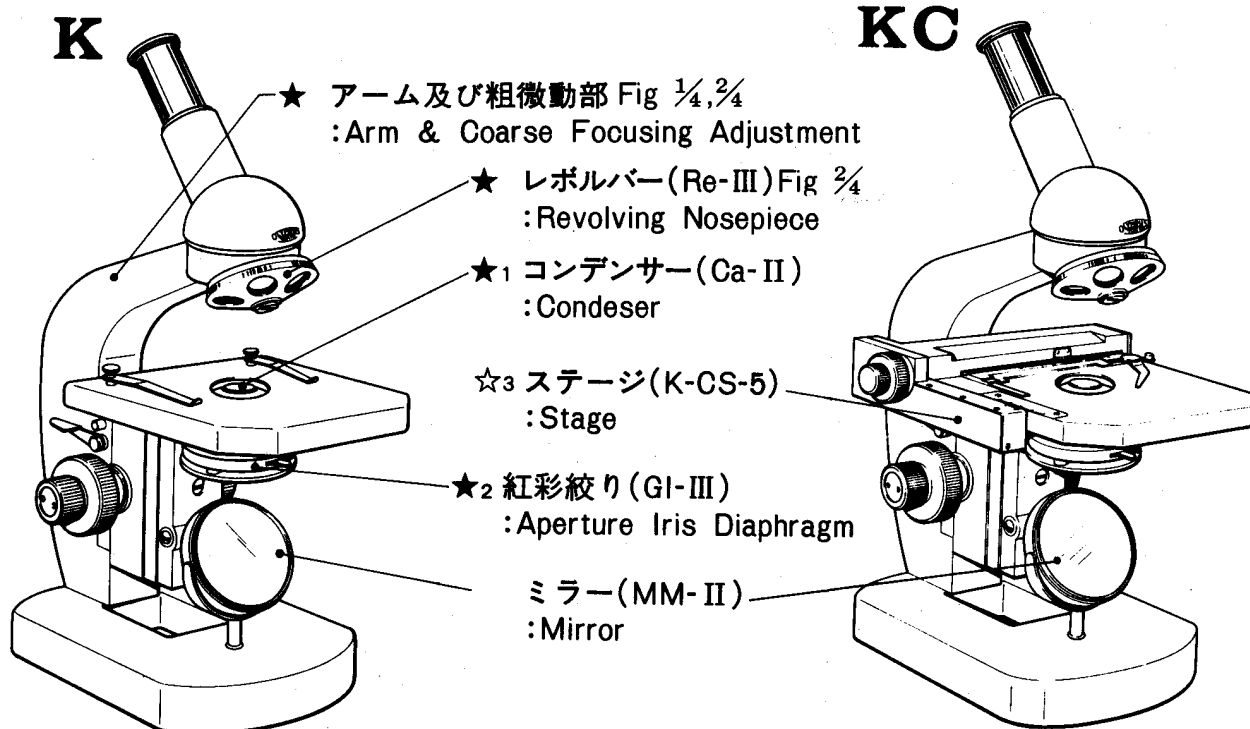


**K**

**KC**

- ★ アーム及び粗微動部 Fig 1/4, 2/4  
: Arm & Coarse Focusing Adjustment
- ★ レボルバー (Re-III) Fig 2/4  
: Revolving Nosepiece
- ★1 コンデンサー (Ca-II)  
: Condenser
- ☆3 ステージ (K-CS-5)  
: Stage
- ★2 紅彩絞り (GI-III)  
: Aperture Iris Diaphragm
- ミラー (MM-II)  
: Mirror



**KHC-Bi**

**KHS-Bi/mono**

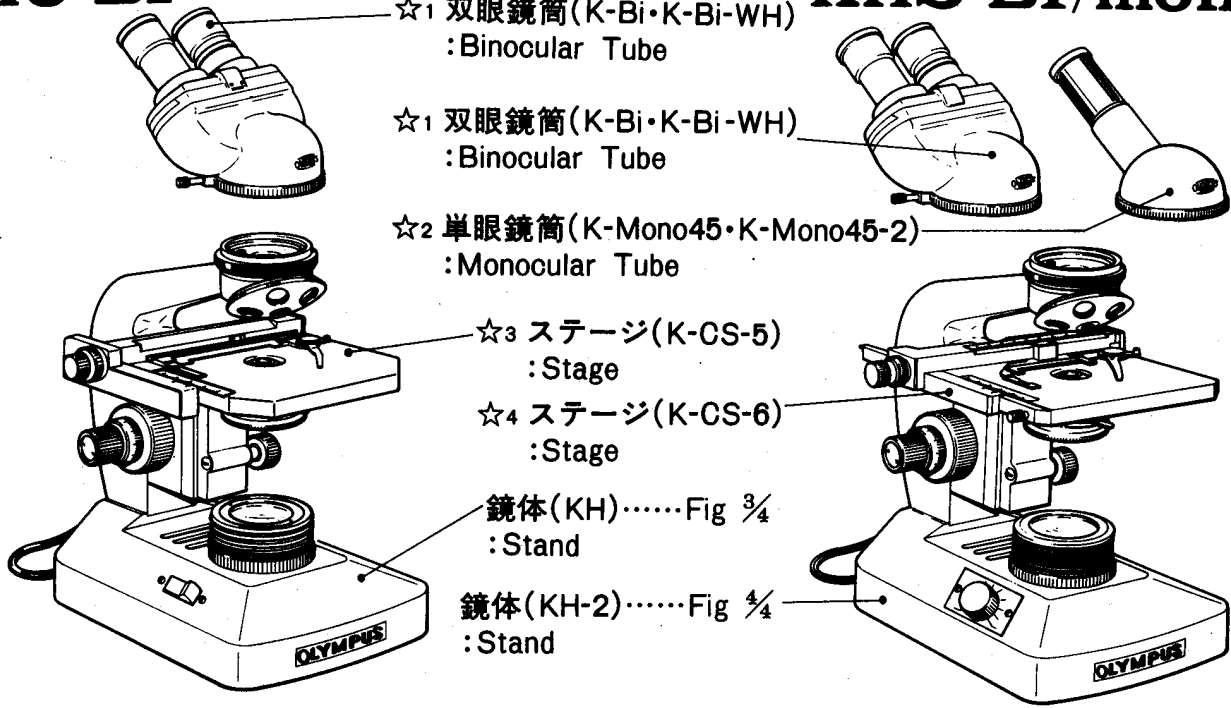

- ☆1 双眼鏡筒 (K-Bi・K-Bi-WH)  
: Binocular Tube
- ☆1 双眼鏡筒 (K-Bi・K-Bi-WH)  
: Binocular Tube
- ☆2 単眼鏡筒 (K-Mono45・K-Mono45-2)  
: Monocular Tube
- ☆3 ステージ (K-CS-5)  
: Stage
- ☆4 ステージ (K-CS-6)  
: Stage
- 鏡体 (KH) ..... Fig 3/4  
: Stand
- 鏡体 (KH-2) ..... Fig 4/4  
: Stand

(注) ★印は、一部の部品を除き、Kシリーズ全機種に共通。: Common parts in K Series with some exceptions.  
 ★1. ★2「12 顕微鏡附属品」を参照。Refer to PARTS LIST No.12.  
 ☆印は、一部の部品を除き、他機種と共通。: Common parts with other microscope models with some exceptions.  
 ☆1「12 顕微鏡附属品」のBi-H45-CH-2を参照。: Refer to Model Bi-H45-CH-2 in PARTS LIST No.1  
 ☆2「12 顕微鏡附属品」のMono-45あるいはMono-45-2を参照。: Refer to Mono-45 or Mono-45-2 in PARTS LIST No.12.  
 ☆3「12 顕微鏡附属品」のCS-5を参照。: Refer to Model CS-5 in PARTS LIST No.12.  
 ☆4「12 顕微鏡附属品」のCS-6を参照。: Refer to Model CS-6 in PARTS LIST No.12.

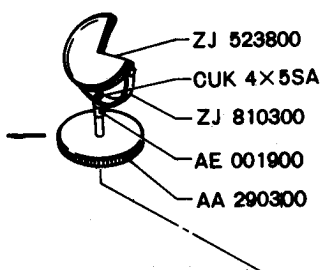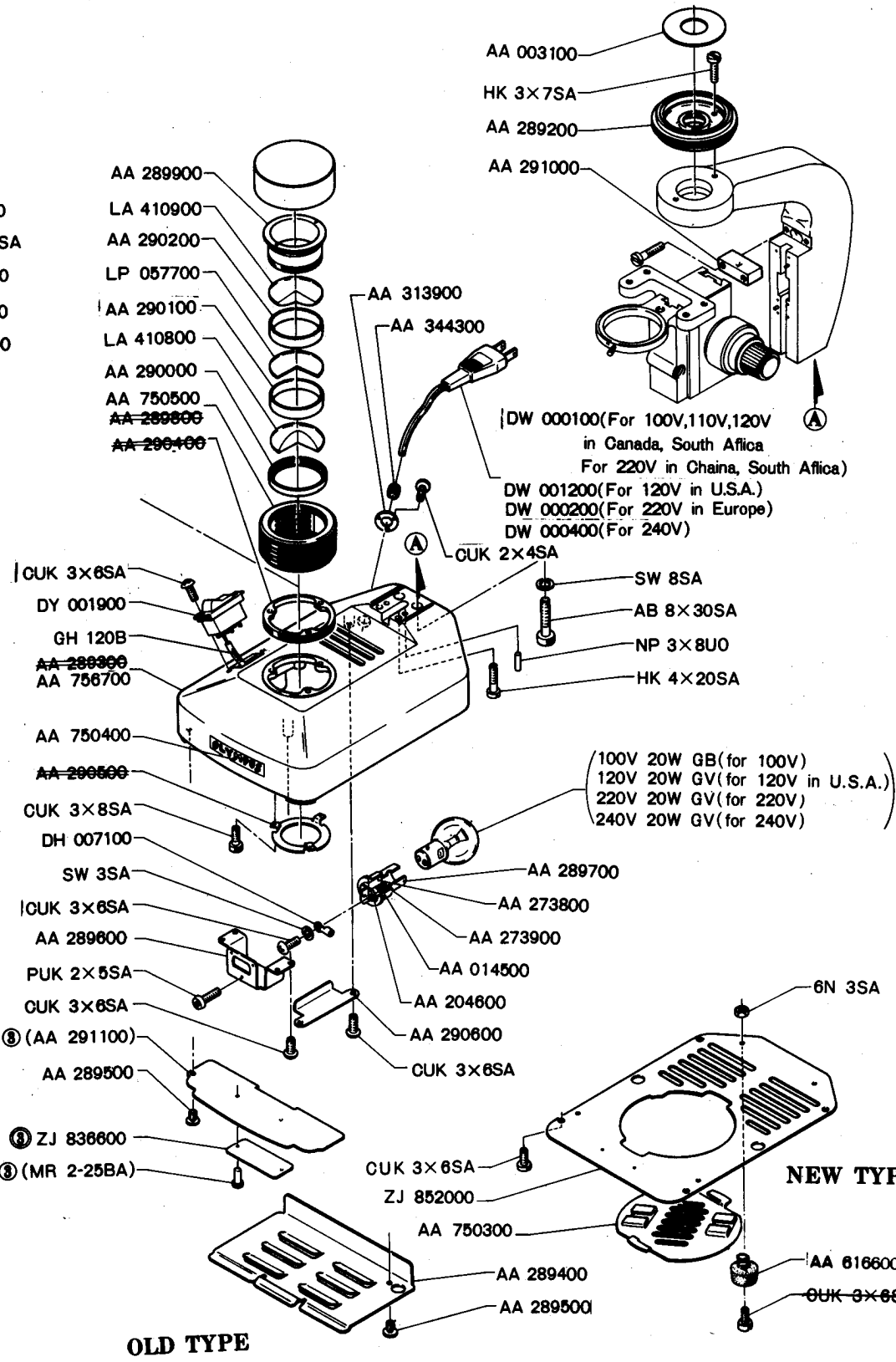
**A****B****C****D****1****MM-II****2****3****4****EXPLODED PARTS DIAGRAM**

MODEL	UNIT	FIG
K Series	KH	3/4

**OLYMPUS OPTICAL CO., LTD. TOKYO, JAPAN**

NOTE: WHEN ORDERING FOR SPARE PARTS, PLEASE CLARIFY A MODEL OR UNIT, PARTS NUMBER AND QUANTITY.

②

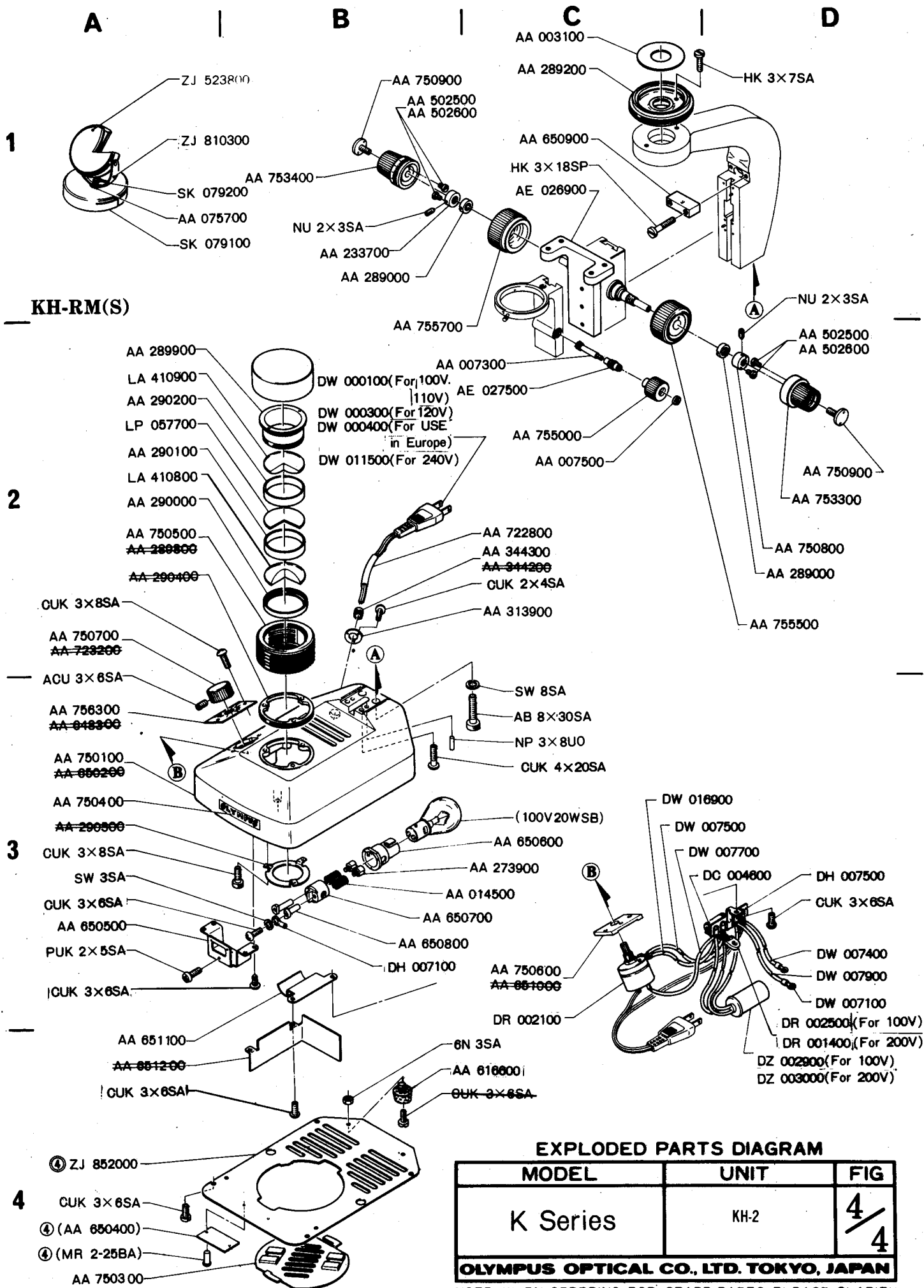

**EXPLODED PARTS DIAGRAM**

MODEL	UNIT	FIG
K Series	KH-2	4/4

**OLYMPUS OPTICAL CO., LTD. TOKYO, JAPAN**

NOTE: WHEN ORDERING FOR SPARE PARTS, PLEASE CLARIFY A MODEL OR UNIT, PARTS NUMBER AND QUANTITY.

結線図 ; CONNECTING DIAGRAM

